

HAMAMATSU

DATA SHEET

Electron Multiplier CCD Digital Camera



C9100-02 (1000 × 1000)

C9100-12 (BT-CCD)




The C9100 cameras bring together all the advantages of the latest Electron Multiplication CCDs with Hamamatsu engineering. High gain, good signal to noise, resolution and speed are combined with a proprietary hermetic vacuum chamber evacuated to 10^{-8} Torr. High vacuum, deep cooling and specially designed electronics combine to reduce camera noise before the gain multiplication process begins.

Using on-chip electron gain multiplication and deep cooling, the signal level can be greatly increased relative to the camera noise; providing a relative read noise of less than one electron at high gain levels. Depending on the model, gain factors up to 800 times or 2000 times are possible while still operating at 30 frames per second at full spatial resolution and 14 bit digitization. Frame rates of 250 frames per second or greater are possible by using the binning and sub-array features of the C9100 cameras.

The C9100 series cameras are recommended for any application requiring, speed, signal to noise, dynamic range and resolution at low light levels.

SPECTRAL RESPONSE CHARACTERISTICS

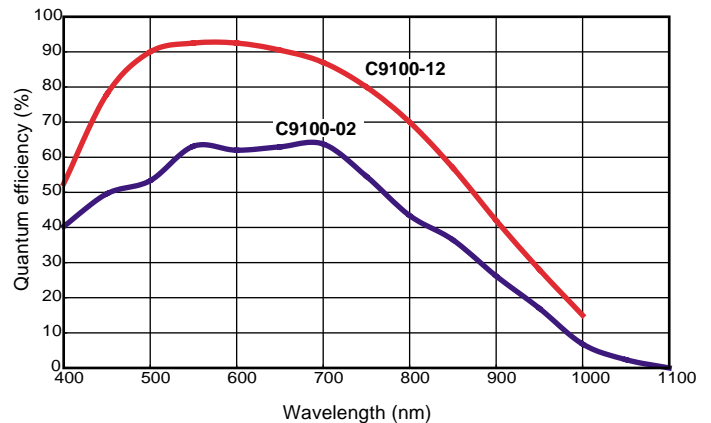

★ This is typical, not guaranteed


FEATURES

- High EM gain
 - 800 times for C9100-02
 - 2000 times for C9100-12
- - 50 °C cooling with hermetic sealed head
- Real time (> 30 Hz) readout with full resolution
 - 35 MHz clock for C9100-02 (1000 × 1000)
 - 11 MHz clock for C9100-12 (512 × 512)

APPLICATIONS

- Real time imaging for low light imaging
- Real time confocal microscopy
- I.I. readout for single molecule fluorescence imaging
- TIRF microscope imaging

SYSTEM CONFIGURATION

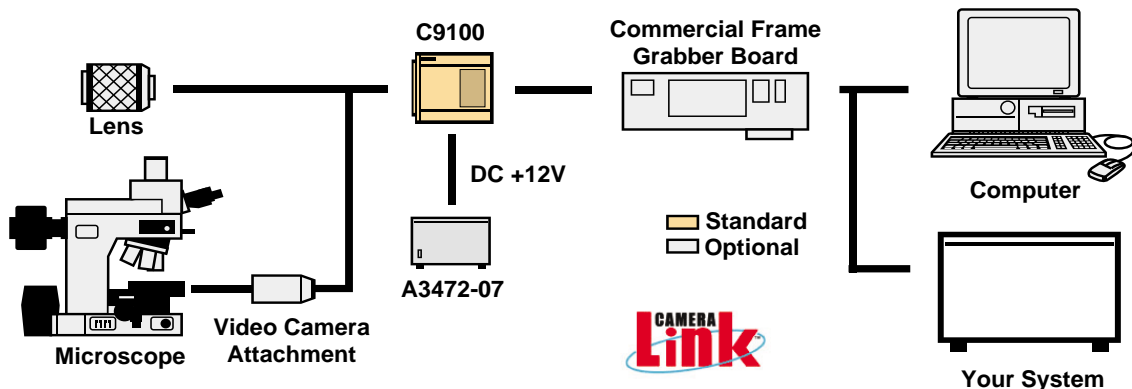

The C9100 camera outputs data is transmitted via a Camera Link interface.

SPECIFICATIONS


Type number	C9100-02	C9100-12		
Camera head type	Hermetic vacuum-sealed air-cooled head			
Circulating water cooler	-			
Mechanical shutter	-			
Imaging device	Frame transfer CCD	Backthinned Frame transfer CCD		
Effective no. of pixels	1000 (H) × 1000 (V)	512 (H) × 512 (V)		
Cell size	8 μm (H) × 8 μm (V)	16 μm (H) × 16 μm (V)		
Effective area	8.0 mm (H) × 8.0 mm (V)	8.192 mm (H) × 8.192 mm (V)		
Pixel clock rate	35 MHz/pixel	11 MHz/pixel		
Frame rate***	1 × 1	30.6 frame/s	35.8 frame/s	
	binning	2 × 2	57.9 frame/s	66.9 frame/s
		4 × 4	104.6 frame/s	118.2 frame/s
		8 × 8	175.1 frame/s	191.6 frame/s
		16 × 16	264.1 frame/s	277.0 frame/s
Readout noise(r.m.s.) typ.	at EM-gain min.	30 electrons	100 electrons	
	at EM-gain max.	< 1 electrons	< 1 electrons	
Full well capacity typ.	70000 electrons	400000 electrons		
Electron multiplier gain max.	× 800*	× 2000*		
Cooling method	Forced-air peltier cooling with hermetic sealing**			
Cooling temperature	- 50 °C			
A/D converter	14 bit			
Output signal / External control	Camera Link			
Exposure time	100 μs to 10 s	27.1 ms to 10 s		
Electronic shutter	Yes	No		
Sub-array***	Yes			
External trigger	Yes			
Offset enhancement	Yes			
Contrast enhancement	No			
Lens mount	C-mount			
Power requirements	DC +12 V			
Power consumption	approx. 60 V·A			
Ambient storage temperature	- 10 °C to + 50 °C			
Ambient operating temperature	0 °C to + 40 °C			
Ambient operating/storage humidity	70 % max. (with no condensation)			

* Even with electron multiplier gain maximum, dark signal is kept low level for low light imaging.

** The hermetic sealed head maintains a high degree of vacuum 10⁻⁸ Torr, without re-evacuation.

*** Frame rate of each binning and sub-array condition	Binning	Effective vertical width (Sub-array)												
		C9100-02						C9100-12						
		1000	512	256	128	64	32	16	512	256	128	64	32	16
Frame rate (frame/s)	1 × 1	30.6	56.7	102.6	172.3	260.9	351.2	424.7	35.8	66.9	118.3	191.9	279.3	361.0
	2 × 2	57.9	102.6	172.3	260.9	351.2	424.7	474.4	66.9	118.3	191.9	279.3	361.0	421.9
	4 × 4	104.6	172.3	260.9	351.2	424.7	474.4	503.8	118.2	191.9	278.6	359.7	421.9	460.8
	8 × 8	175.1	260.9	351.2	424.7	474.4	503.8	520.0	191.6	278.6	359.7	420.2	460.8	483.1
	16 × 16	264.1	351.2	424.7	474.4	53.8	520.0	528.4	277.0	357.1	418.4	456.6	478.5	490.2

DIMENSIONAL OUTLINES (Unit: mm)

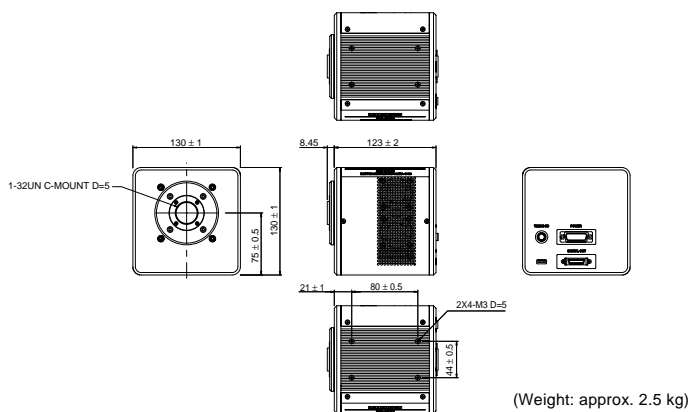

OPTIONAL

- **Power supply unit: A3472-07**
Line voltage: 100 to 240V AC input
Output voltage: DC +12V
Dimension: 182 mm (W) × 240 mm (D) × 60 mm(H)
- **Power cable: A9189-05**



★ Windows is registered trademark of Microsoft Corporation in the U.S.A.

★ Product and software package names noted in this documentation are trademarks or registered trademarks of their respective manufacturers.

- Subject to local technical requirements and regulations, availability of products included in this promotional material may vary. Please consult with our sales office.
 - Information furnished by HAMAMATSU is believed to be reliable. However, no responsibility is assumed for possible inaccuracies or omissions.
- Specifications and external appearance are subject to change without notice.

© 2003 Hamamatsu Photonics K.K.

HAMAMATSU

Homepage Address <http://www.hamamatsu.com>

HAMAMATSU PHOTONICS K.K., Systems Division

812 Joko-cho, Hamamatsu City, 431-3196, Japan, Telephone: (81)53-431-0124, Fax: (81)53-435-1574, E-mail: export@sys.hpk.co.jp

U.S.A. and Canada: Hamamatsu Photonic Systems: 360 Foothill Road, Bridgewater, N.J. 08807-0910, U.S.A., Telephone: (1)908-231-1116, Fax: (1)908-231-0852, E-mail: usa@hamamatsu.com

Germany: Hamamatsu Photonics Deutschland GmbH: Arzbergerstr. 10, D-82211 Herrsching am Ammersee, Germany, Telephone: (49)8152-375-0, Fax: (49)8152-2658, E-mail: info@hamamatsu.de

France: Hamamatsu Photonics France S.A.R.L.: 8, Rue du Saule Trapu, Parc du Moulin de Massy, 91882 Massy Cedex, France, Telephone: (33)1 69 53 71 00, Fax: (33)1 69 53 71 10, E-mail: info@hamamatsu.fr

United Kingdom: Hamamatsu Photonics UK Limited: 2 Howard Court, 10Tewin Road, Welwyn Garden City, Hertfordshire, AL7 1BW, U.K., Telephone: (44) 1707-294888, Fax: (44) 1707-325777, E-mail: info@hamamatsu.co.uk

North Europe: Hamamatsu Photonics Norden AB: Smidesvägen 12, SE-171-41 Solna, Sweden, Telephone: (46)8-509-031-00, Fax: (46)8-509-031-01, E-mail: info@hamamatsu.se

Italy: Hamamatsu Photonics Italia S.R.L.: Strada della Mois, 1/E 20020 Arese (Milano), Italy, Telephone: (39)02-935 81 733, Fax: (39)02-935 81 741, E-mail: info@hamamatsu.it

Cat. No. SICS1123E01
NOV/2003 HPK
Created in Japan (PDF)